

GET YOUR SPILLS UNDER CONTROL WITH OUR GREEN-Z SPILL CONTROL SOLIDIFIERS



Solidify Chemical Spills, Blood, and Bodily Fluid Spills with Ease.

Green-Z® Spill Control Solidifier | Chlorine-Free Formula

At times, workers may be unintentionally exposed to blood and bodily fluid spills along with other harmful substances such as toxic chemicals, biological hazards and chemotherapy agents. To help reduce those exposure risks, rely on Green-Z® for improved safety in your spill cleanup process.



Fast Acting



Chlorine Free



Splash Free

Features

- Contains fast-acting, superabsorbent polymers
- Chlorine-free (safe to use on chemotherapy, glutaraldehyde, and mercury spills)
- Turns spills into semi-solid (gel) mass
- Cost-effective cleanup solution

Directions for Use

1. Wear appropriate protective clothing.
2. Sprinkle Green-Z powder to surround the spill and then evenly distribute over it. Allow sufficient time for solidification to occur.
3. Remove solidified material with scoop and scraper (or similar tool) and place in an appropriate container.
4. Clean and disinfect contaminated areas after use and wash thoroughly.

 **PharmaSystems®**

PharmaSystems.com
151 Telsons Road
Markham Ontario
Canada L3R 1E7

Item No.: 42012

Size: 15 oz.

Case Qty: 12

Expiration Dating

Green-Z® is shelf stable and does not require an expiration date. In order for the product to remain shelf stable, it must be kept in a cool, dry place and in its original packaging. If contents are hard or compact, simply shake the product to loosen, before use.

